

PRODUCT DATA DIPSLIDE DISPOSAL

Product Uses

For the monitoring of surfaces and liquids for microbiological contamination.

Product Appearance

Plastic tube containing a paddle coated on both sides with a nutritive gel.

Product Storage

Cool dry place not refrigerated.

Product Fire Risk

If incinerated, plastic components will burn and may produce toxic fumes.

Product Corrosiveness

None.

Product Irritation

None known

Product Precautions

Used dipslides may contain viable micro-organisms. Avoid opening the tubes, DO NOT ingest or touch the contaminated gel.

Product Disposal

Sterilise before disposal. Incinerate, autoclave at 121⁰C for 15 minutes or immerse overnight in 25% solution of household bleach or 5% solution of phenolic disinfectant. Sterilised dipslides may be disposed of in the normal way.